

The genie in the bottle

Education in Chemistry

September 2018

rsc.li/2PyWj3w



Technician notes

Kit

- Approx. 50 cm³ of ethanol (importantly, not IDA/IMS), (highly flammable liquid and vapour, harmful) (*Hazard 40A, recipe 002*)
- Boric acid (may damage fertility, may damage the unborn child) (*Hazard 14A*)
- 2.5 – 3.0 g of the following salts:
 - lithium chloride (*Hazard 47B*)
 - sodium chloride (*Hazard 47B*)
 - potassium chloride (*Hazard 47B*)
 - anhydrous calcium chloride (irritant) (*Hazard 19A*)
 - hydrated strontium chloride (*Hazard 19A*)
 - hydrated barium chloride (harmful, toxic) (*Hazard 10A*)
 - hydrated copper(II) chloride (harmful, irritant, very toxic to aquatic life) (*Hazard 27A*)
- 80 mm borosilicate crystallising dishes or 250 cm³ borosilicate beakers
- Heat-resistant mats
- Wooden splints
- Metre rule

Preparation

Work in a well-ventilated room and wear eye protection. Place the beakers / basins on the heat-resistant mats and spread 2.5 – 3.0 g of each salt around the base of the container. Dampen the solid with approximately 0.5 cm³ of water and add approximately 6 cm³ ethanol over each. Stopper the ethanol bottle and remove it to at least 2 m away from the demonstration.

In front of the class

Position the audience 3 m away from the demonstration with eye protection. Ask for a volunteer to switch off the lights if the switch is not nearby. Ignite the solvent in each container with a splint on a metre rule. Each beaker will burn with a characteristic colour. Allow the flames to burn out – do not attempt to add more solvent until at least 15 minutes have passed since the last flame extinguished.

Safety and disposal

Boric acid is a teratogen – you may not wish to use it.

Do not be tempted to pour more alcohol onto the flame or hot glassware in order to extend the demonstration. Use only borosilicate glass (avoid watchglasses, which are usually made from soda glass and may crack).

Dilute the contents of the copper(II) chloride beaker to 1 L and pour down the sink. The other salts will be unchanged at the end of the demonstration and can be retrieved and recycled for future demonstrations.